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Table1 Antigens and antibodies used

Abbreviation of analyte antibody	Name of Antibody/Antigen	Specification	Supplier
MAb1	MAB287	Mouse IgG1	R&D Systems, Inc.
MAb2	MAB1072	Mouse MAb	Chemicon (Millipore Corp)
MAb3	BM802	Mouse IgG1	Acris Ab GmbH
Poly1	AB-286-NA	Total Rabbit IgG	R&D Systems, Inc.
Poly2	LS-C-11323	Rabbit Polyclonal (IgG)	Life Span Bioscience
EPO	ESPO	Recombinant Human Erythropoietin- α	Kyowa Hakko Kirin Pharma.
	EPO- α		Prospec-Tany Techno gene Ltd.

Table2 Assay Condition

Assay format	Drug	Detection antibody
Direct coating method	ESPO	Use ①* for MAb1 to 3 Use ②** for Poly1 and 2
Indirect coating method	Biotin labeled EPO- α	Use ①* for MAb1 to 3
Bridging assay	1. Biotin labeled EPO- α 2. SULFO-TAG labeled EPO- α	Not used

① * : SULFO-TAG labeled goat anti-mouse IgG antibody

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Table3 Detection range

Abbreviation of analyte antibody	Direct coating method	Indirect coating method	Bridging assay
MAb1	78.1~5,000 ng/mL	0.012~500 ng/mL	4.88~2,500 ng/mL
MAb2	Not detected	0.049~200 ng/mL	4.88~2,500 ng/mL
MAb3	Not detected	0.2~200 ng/mL	4.88~2,500 ng/mL
Poly1	39.1~5,000 ng/mL		39.1~2,500 ng/mL
Poly2	39.1~5,000 ng/mL		39.1~2,500 ng/mL

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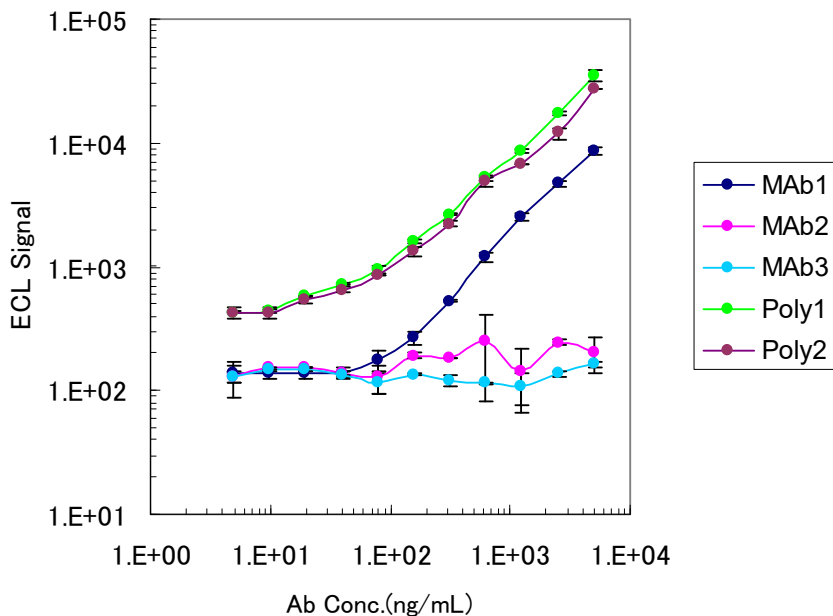
Fig. 1 Calibration curve of direct coating method

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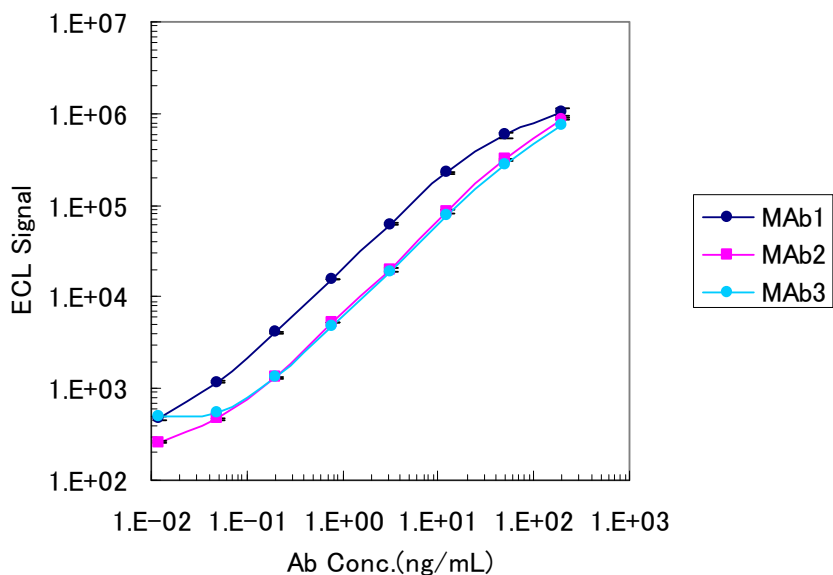
Fig.2 Calibration curve of indirect coating method

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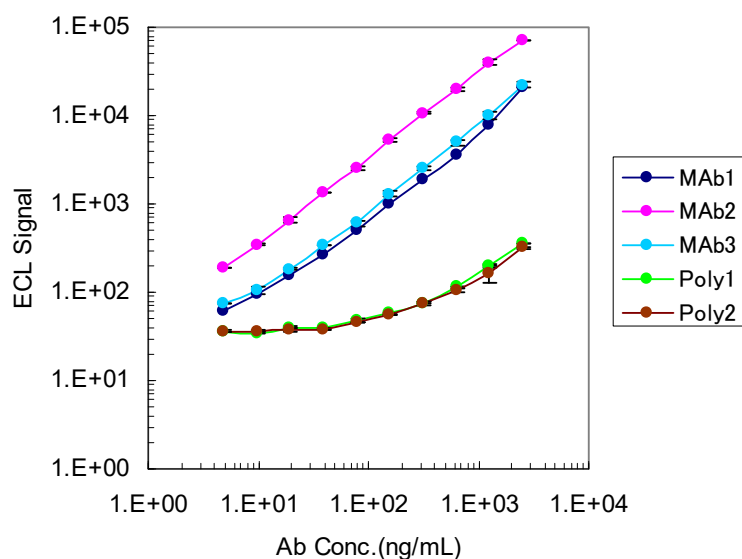
Fig.3 Calibration curve of bridging assay

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- 1) Anthony R. Mire-Sluis, Yu Chen Barrett, Viswanath Devanarayan, Eugen Koren, Hank Liu, Mauricio Maia, Thomas Parish, George Scott, Gopi Shankar, Elizabeth Shores, Steven J. Swanson, Gary Taniguchi, Daniel Wierda, Linda A. Zuckerman, *J. Immunol. Methods*, 289, 1(2004).
- 2) Guideline on Immunogenicity assessment of therapeutic proteins, European Medicines Agency, Committee for Medicinal Products for Human Use (CHMP), May 2017
- 3) Meso Scale Discovery, Bridging Immunogenicity Assays Guidelines for Assay Development
<<https://www.mesoscale.com/~media/files/technical%20notes/bridging%20immunogenicity%20assay.pdf?la=ja-jp>>(accessed 2025.7.2)

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